

WCSB, Wageningen University	IBISBA-SOP-WU13
	Version 1.0

EPP - Standard Operating Procedure

(only for selected experiments intended to transfer results from one lab to the other)

Title: Colony PCR of *S. cerevisiae*, *P. putida*, *E. coli*

<u>distribution list</u>			
changes to prior version:			
	name	signature	date
Experimenter 1	Rita Volkers		14/3/2019

Instruction

Colony PCR of *S. cerevisiae*, *P. putida*, *E. coli*

1. Introduction / Purpose

In this SOP, colony PCR of *S. cerevisiae*, *P. putida*, and *E. coli* is described. There are alternative methods available for both bacterial strains that are not described here.

Keywords: Colony – PCR – *S. cerevisiae* – *P. putida* – *E. coli*

2. Equipment and chemicals

2.1. Equipment

- PCR machine

2.2. Chemicals

- Phire Plant Direct PCR Kit (Thermo Fisher Scientific)
- Phire Green Hot Start II DNA Polymerase (Thermo Fisher Scientific)
- MQ water

2.3 Other materials

- Primers (forward and reverse)
- The colonies have to be analysed by PCR

3. Procedures

The template DNA in a colony PCR reaction is prepared as follows:

- Put 30 µl of Dilution Buffer (from Phire Plant Direct PCR Kit) in a 1.5 ml eppendorf tube
- Transfer a colony from an agar plate to the tube, resuspend
- Incubate 2 minutes @ 70°C in a heat block
- Use 1 µl of this mixture as template

Prepare PCR reaction mixture with Phire Green Hot Start II DNA Polymerase kit:

- MQ water: enough to make a total volume of 20µl
- 4 µl 5x Phire Green Reaction Buffer
- 0.4 µl 10 mM dNTP's

- 1 μ l Primer forward(stock of 10 pmol/ μ l)
- 1 μ l Primer reverse (stock of 10 pmol/ μ l)
- 0.4 μ l Phire Hot Start II DNA Polymerase
- 1 μ l DNA template

Volume can be scaled up (to 50 μ l) or down (to 15 μ l), change all volumes accordingly.

Cycling protocol:

Cycling Step	Temperature	Time	Cycles
Initial denaturation	98°C	30 s	1
Denaturation Annealing Extension	98°C	5 s	25-35
	X°C	5 s	
	72°C	10-15 s/kb	
Final extension	72°C	1 min	1
Hold	16°C	hold	1

Annealing temperature depends on primers.

5. Remarks / troubleshooting

Because of the use of 5x Phire Green Reaction Buffer, addition of loading dye is not necessary before running an agarose gel.

Troubleshooting: see manuals of kits on website of Thermo Fisher Scientific.

6. Biosafety

No biosafety issues are associated with this protocol.

7. Acknowledgements



IBISBA (grant agreement No 730976) and EmPowerPutida (grant agreement No 635536) projects have received funding from the European Union's Horizon 2020 research and innovation programme.